

# SOP for Colour Measurement in Fresh Yam (*Dioscorea* Sp.) and Fresh Cassava (*Manihot esculenta*) using Chromameter

High-Throughput Phenotyping Protocols (HTPP), WP3

Lusaka, Zambia, July 2020

Emmanuel ALAMU, International Institute of Tropical Agriculture (IITA), Lusaka, Zambia

Michael ADESOKAN, IITA, Ibadan, Nigeria

Karima MEGHAR, Centre de coopération Internationale en Recherche Agronomique pour le Développement (CIRAD), Montpellier, France (Validator)

Fabrice DAVRIEUX, CIRAD, Saint-Pierre, France-Réunion (Validator)



This report has been written in the framework of RTBfoods project.


To be cited as:

**Emmanuel ALAMU, Michael ADESOKAN, Karima MEGHAR, Fabrice DAVRIEUX** (2020). *Standard Operating Protocol for Colour Measurement in Fresh Yam (*Dioscorea* Sp.) and Fresh Cassava (*Manihot esculenta*) using Chromameter. High-Throughput Phenotyping Protocols (HTPP), WP3*. Lusaka, Zambia: RTBfoods Laboratory Standard Operating Procedure, 10 p. <https://doi.org/10.18167/agritrop/00664>

Ethics: The activities, which led to the production of this manual, were assessed and approved by the CIRAD Ethics Committee (H2020 ethics self-assessment procedure). When relevant, samples were prepared according to good hygiene and manufacturing practices. When external participants were involved in an activity, they were priorly informed about the objective of the activity and explained that their participation was entirely voluntary, that they could stop the interview at any point and that their responses would be anonymous and securely stored by the research team for research purposes. Written consent (signature) was systematically sought from sensory panelists and from consumers participating in activities.

Acknowledgments: This work was supported by the RTBfoods project <https://rtbfoods.cirad.fr>, through a grant OPP1178942: Breeding RTB products for end user preferences (RTBfoods), to the French Agricultural Research Centre for International Development (CIRAD), Montpellier, France, by the Bill & Melinda Gates Foundation (BMGF).

Image cover page © LAJOUS P. for RTBfoods.

<b>RTBfoods</b> <b>WP3: High-throughput phenotyping protocols (HTPP)</b>		
<b>SOP: Colour Measurement in Fresh Yam (Dioscorea Sp.) and Fresh cassava (Manihot esculenta) using Chromameter</b>		
<b>Date: 08/07/2020</b>		<b>Release: 2</b>
<p>Written by: Emmanuel Oladeji ALAMU &amp; Michael ADESOKAN</p> <p>For information on this SOP please contact:</p> <ul style="list-style-type: none"> <li>• Busie MAZIYA-DIXON, <a href="mailto:b.maziya-dixon@cgiar.org">b.maziya-dixon@cgiar.org</a></li> <li>• Emmanuel Oladeji ALAMU, <a href="mailto:oalamu@cgiar.org">oalamu@cgiar.org</a></li> <li>• Michael ADESOKAN, <a href="mailto:m.adesokan@cgiar.org">m.adesokan@cgiar.org</a></li> </ul>		
<p align="center"><b>This document has been approved by:</b></p>		
Partner	Name of the person who approved	Date
CIRAD	Karima MEGHAR	08/07/2020
CIRAD	Fabrice DAVRIEUX	09/07/2020

# CONTENTS

---

## Table of Contents

1	Scope and Application .....	5
2	References .....	5
3	Principle.....	5
4	Reagents .....	6
5	Apparatus .....	6
6	Procedure .....	8
7	Expression of Results .....	9
8	Critical Points or Note on the Procedure .....	9
9	Revision Record .....	9

**RTBfoods- WP3: High-throughput phenotyping protocols (HTPP)**  
**SOP: Colour Measurement in Fresh Yam (*Dioscorea* Sp.) and Fresh cassava (*Manihot esculenta*) using Chromameter**

**Date: 08/07/2020**

**Release: 2**

## 1 SCOPE AND APPLICATION

Colour is a characteristic of light, which is measurable in terms of intensity and wavelength. The primary pigments imparting colour quality are the fat-soluble chlorophylls (green), carotenoids (yellow, orange, and red), water-soluble anthocyanins (red, blue), flavonoids (yellow) and betalains (red) (Barrett et al. 2010). The measurement of Colour could be determined by human vision, which could be subjective. Thus, colour measurement using the instrument is recommended. Colour can be correlated with other quality attributes such as sensory, nutritional, and visual or non-visual defects. Enzymatic browning of yam caused by the oxidation of phenols by polyphenol oxidases and peroxidases is a common phenomenon associated with yam discolouration during injury or processing at low temperature. The extent of this discolouration is determined by the  $L^*$ ,  $a^*$  and  $b^*$  values of measurement using the Chromameter.

## 2 REFERENCES

Barrett, D. M., Beaulieu, J. C., & Shewfelt, R. (2010). Colour, flavour, texture, and nutritional quality of fresh-cut fruits and vegetables: desirable levels, instrumental and sensory measurement, and the effects of processing. *Critical Reviews in Food Science and Nutrition*, 50(5), 369–389.

Granato, D., & Masson, M. L. (2010). Instrumental Colour and sensory acceptance of soy-based emulsions: a response surface approach. *Ciência e Tecnologia de Alimentos*, 30(4), 1090–1096.

Hutchings, J. B. (1999) *Food Colour and Appearance* (London: Chapman and Hall Publisher)

Pankaj BP and Umezuruile L.O. 2013. Colour Measurement and Analysis in Fresh and Processed Foods: A Review. *Food and Bioprocess Technology*. vol: 6 pages 36-60. DOI: 10.1007/s11947-012-0867-9.

## 3 PRINCIPLE

The CIELAB colour scale is approximately uniform, and the differences between points plotted in the colour space correspond to visual differences between the colours plotted. The CIELAB colour space is organized in a cube form. The  $L^*$  axis runs from top to bottom. The maximum for  $L^*$  is 100, which represents a perfect reflecting diffuser, the minimum for  $L^*$  is red, Negative  $a^*$  is green, Positive  $b^*$  is yellow and Negative  $b^*$  is blue (Fig 1). However, the HunterLab  $L^*$ ,  $a^*$ ,  $b^*$  values are directly read using portable Minolta CR-400 Chroma Meter.  $L^*$  is an approximate measurement of luminosity, which is the property according to which each Colour can be considered as equivalent to a member of the greyscale, between black and white (Granato and Masson 2010).

**RTBfoods- WP3: High-throughput phenotyping protocols (HTPP)**  
**SOP: Colour Measurement in Fresh Yam (Dioscorea Sp.) and Fresh cassava (Manihot esculenta) using Chromameter**

**Date: 08/07/2020**

**Release: 2**

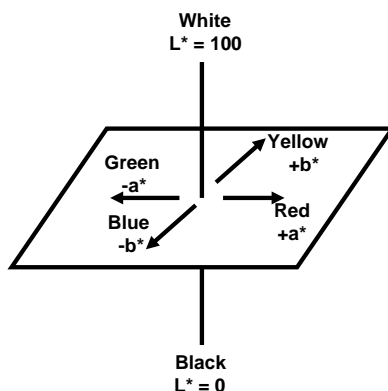




Fig 1. CIELAB colour space

## 4 REAGENTS

No reagent is required

## 5 APPARATUS

Material	Image
<b>Stainless steel knife.</b>  The stainless-steel knife is carbon steel with added chromium to resist corrosion and other elements which increase performance levels.	
<b>Sample trays</b>  Disposable antistatic square weighing trays. Dimension 124 x 22 mm	

**RTBfoods- WP3: High-throughput phenotyping protocols (HTPP)**  
**SOP: Colour Measurement in Fresh Yam (*Dioscorea* Sp.) and Fresh cassava (*Manihot esculenta*) using Chromameter**

**Date: 08/07/2020**

**Release: 2**

Material	Image
<p><b>Mixing Laboratory blender</b></p> <p>The blender is with stainless cup and steel blades</p>	
<p><b>Petri-dishes</b></p> <p><b>Small size 60mm (diameter) x 15mm (height)</b></p> <p>Petri dishes are shallow cylindrical containers with fitted lids and are typically made of borosilicate glass or transparent plastics (usually polystyrene or polycarbonate).</p>	
<p><b>Chroma meter</b></p> <p>(Minolta CR- 400 Chroma Meter)</p>	



**RTBfoods- WP3: High-throughput phenotyping protocols (HTPP)**  
**SOP: Colour Measurement in Fresh Yam (Dioscorea Sp.) and Fresh cassava (Manihot esculenta) using Chromameter**

**Date: 08/07/2020**

**Release: 2**

## 6 PROCEDURE

The Colour of fresh yam samples was measured using Minolta CR- 400 Chroma Meter (Minolta Corp., Osaka, Japan) using L\*a\*b\* systems. It was calibrated against a white standard before taking readings. Colour measurement involves the following steps;

- a) Sampling and sample preparation
- b) Homogenizing and blending of sample
- c) Chromameter reading

The harvest plus sampling protocol was adopted where the fresh yam tuber or fresh cassava root was washed, air-dried, and peeled. A stainless-steel knife was used to cut the tuber/roots into four portions, longitudinally from the proximal to the distal end. Two opposite parts of the four portions were selected and cut into reduced sizes. Chopped pieces were quartered, and two opposite portions were mixed manually to obtain representative sample lots and then transferred into an electric blender (Figure 1). Blended samples were collected in clean, transparent Petri dishes in duplicate and ready for measurement with a Chromameter, as shown in Figure 1.

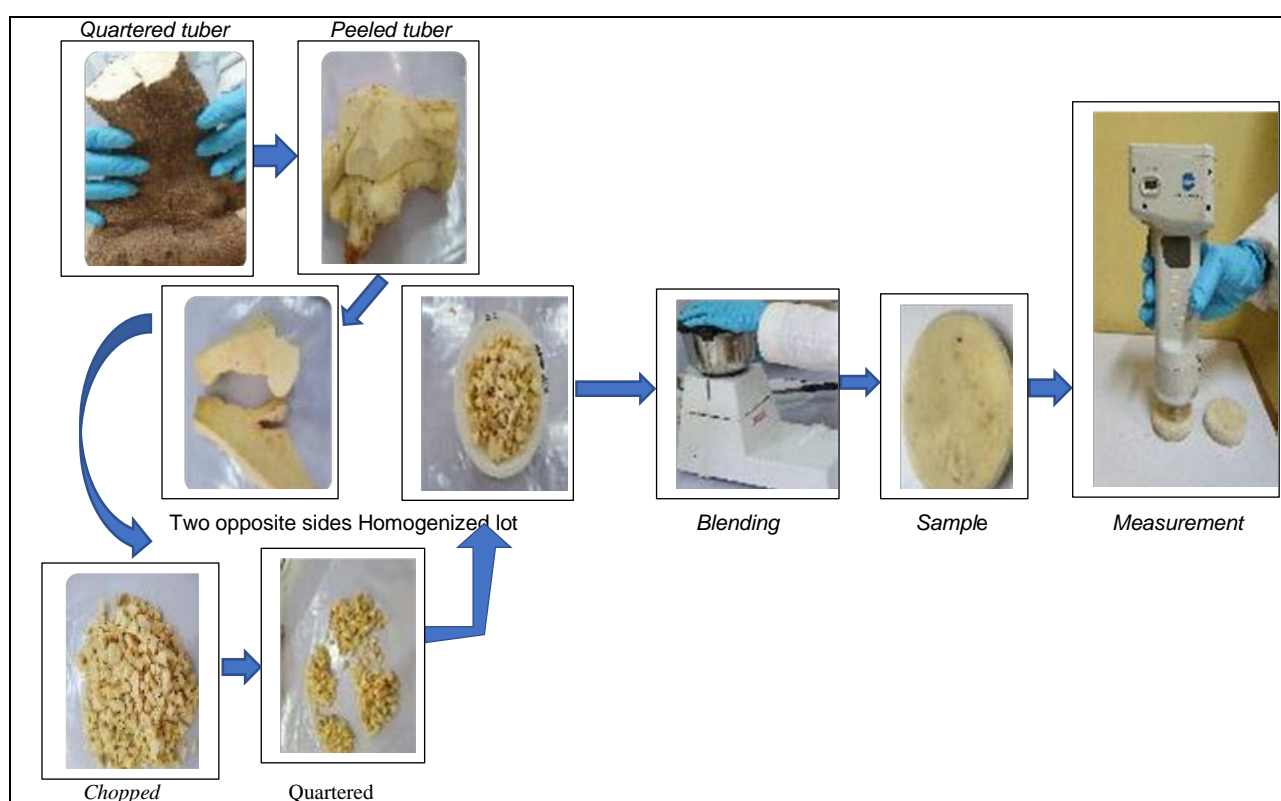


Figure 1: Workflow of color measurements using a Chromameter



**RTBfoods- WP3: High-throughput phenotyping protocols (HTPP)**  
**SOP: Colour Measurement in Fresh Yam (Dioscorea Sp.) and Fresh cassava (Manihot esculenta) using Chromameter**

**Date: 08/07/2020**

**Release: 2**

## 7 EXPRESSION OF RESULTS

**Total colour difference:**

$$\Delta E^2 = L^{2*} + a^{2*} + b^{2*}$$

**Brown Index:** ( Pankaj and Umezurule, 2013)

$$BI = 100 * \frac{(X - 0.31)}{0.17}$$

$$\text{Where } X = \frac{(a^* + 1.75L^*)}{(5.645L^* + a^* - 3.012b^*)}$$

## 8 CRITICAL POINTS OR NOTE ON THE PROCEDURE

- Minolta CR- 400 Chroma Meter must be calibrated against white standard before taking sample readings
- The repeatability test must be carried out by taking at least ten readings on the blended samples. The repeatability uses replicate measurements to determine the standard deviation of the replicates. Chroma values repeatability is an indication of the stability and performance of the Chromameter. The standard deviation (SD) and % Coefficient of variation ( %CV) of the absorbances for ten average spectra are calculated for each wavelength. The SD for L\*, a\* and b\* values of 10 replicates readings were 0.93, 0.25 and 0.63 respectively, while %CV was 1.15, 1.17 and 2.87 respectively. The SD = -1 ≥ 0 ≤ 1 is acceptable. The lower the SD and the % CV ( ≤ 5) the better stability and performance of the Chromameter.
- The sample preparation and reading must be completed within 5 minutes

## 9 REVISION RECORD

Date	Responsible person	Description of change
11/06/2020	Meghar Karima	Reviewing and correction
29/06/2020	Alamu Emmanuel(IITA)	Reviewing, editing and adding input
08/07/2020	Meghar Karima	Reviewing and validation
09/07/2020	Fabrice Davrieux	Validation



**Institution:** Cirad – UMR QualiSud

**Address:** C/O Cathy Méjean, TA-B95/15 - 73 rue Jean-François Breton - 34398  
MONTPELLIER Cedex 5 - France

**Contact Tel:** +33 4 67 61 44 31

**Contact Email:** [rtbfoodspmu@cirad.fr](mailto:rtbfoodspmu@cirad.fr)